Equipment Training Sheet - <u>LECO MSX205M2 Benchtop Sectioning Machine (Cut-off Saw)</u>

Name:		Position:
Departme	ent:	Phone:
		Director:
Ct - CC T	D	
	Trainee Initials	
		 Anyone needing access to MSE equipment needs to contact MSE tech staff: For initial training For subsequent usage (will receive equipment card) No secondary training of the equipment
		 Review of hazards, controls, personal protective equipment: Always wear safety glasses at all times and make sure others are wearing them as well Saw blade spins at several thousand rpm's There are safety interlocks to prevent turning blade on without cover being down Uses 220Volt power
		 Review of standard operating procedure: http://www.mse.engin.umich.edu/internal/procedures/leco-cm-15-cutoff-machine The cutoff saw blades are only available for those associated with the MSE department or classes; otherwise supply your own. Make sure coolant is at a satisfactory level above baffles. Remember to thoroughly clean surfaces with hose provided Lubricate stage, blade and clamp when finished with WD40 Make sure to use appropriate blade Depending on the hardness of material a different blade needs to be used In practice this means that for hard material a "soft" blade is used and gets reduced down faster
		Available supplies to help in the procedure: • Safety glasses • WD-40
O .	ŕ	ate: