


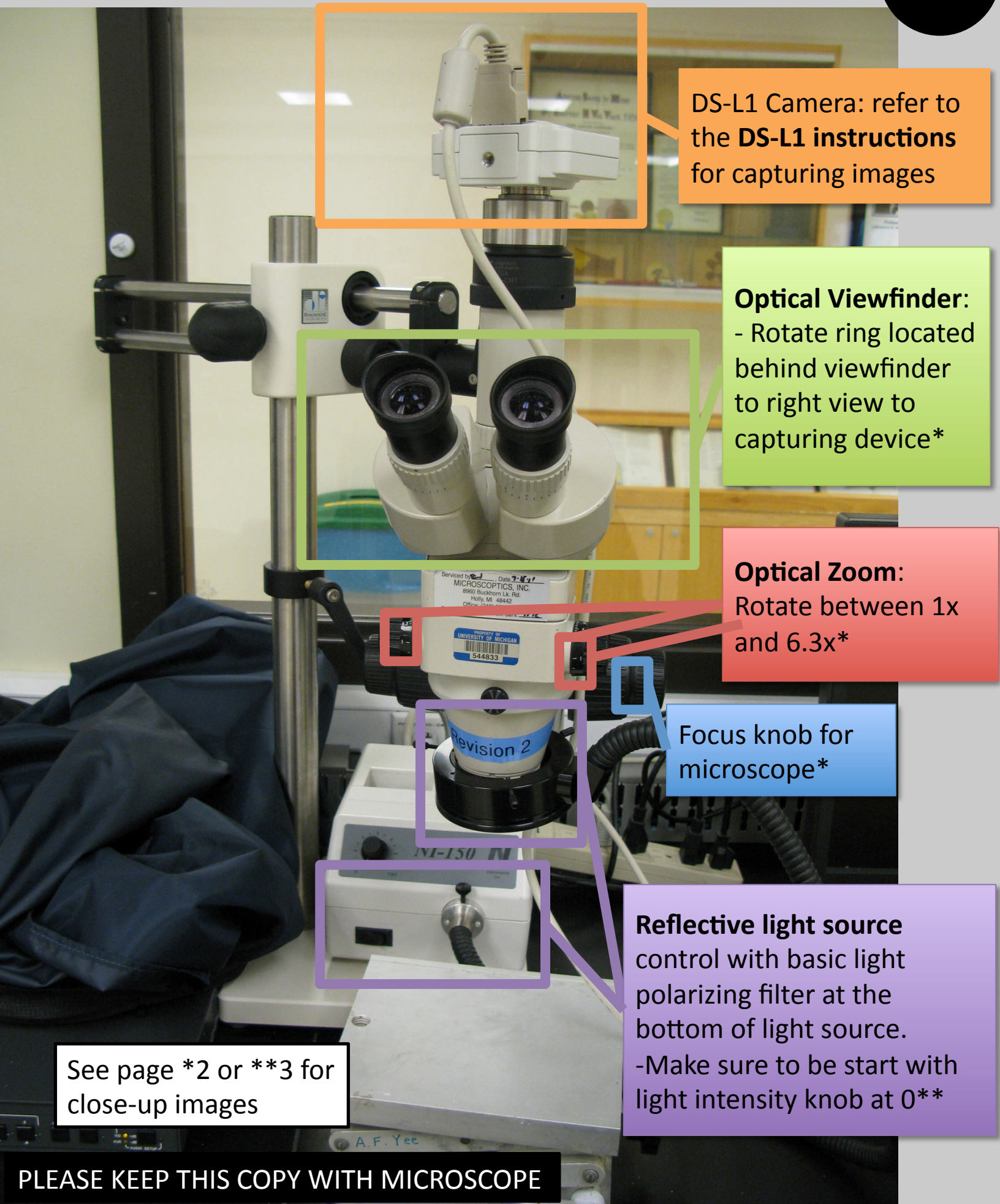

DS-L1 Camera: refer to the **DS-L1 instructions** for capturing images


Optical Viewfinder:
- Rotate ring located behind viewfinder to right view to capturing device*

Optical Zoom:
Rotate between 1x and 6.3x*

Focus knob for microscope*


Reflective light source control with basic light polarizing filter at the bottom of light source.
-Make sure to be start with light intensity knob at 0**

See page *2 or **3 for close-up images

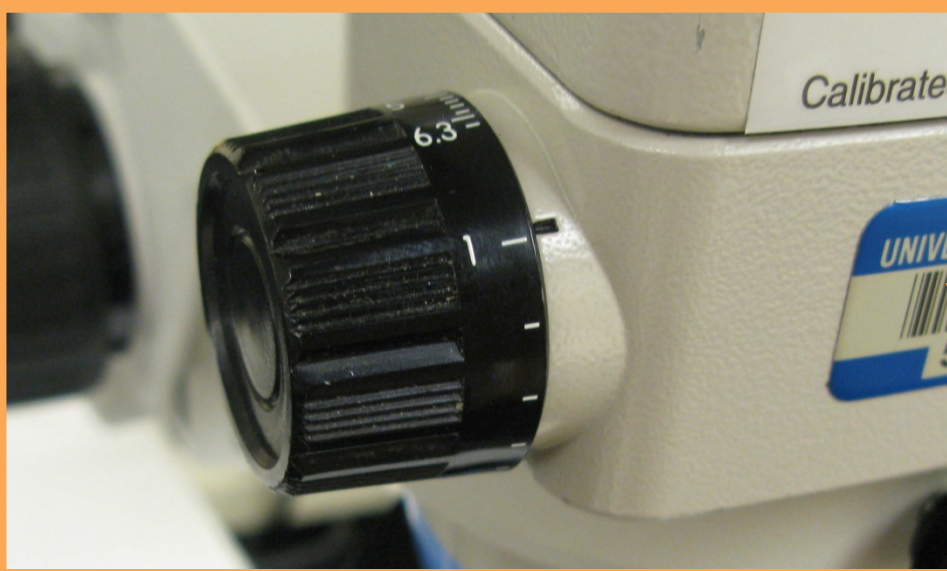


Optical zoom knob with 1x to 6.3x magnification range

Rotate clockwise to transmit light from right eyepiece to capturing device

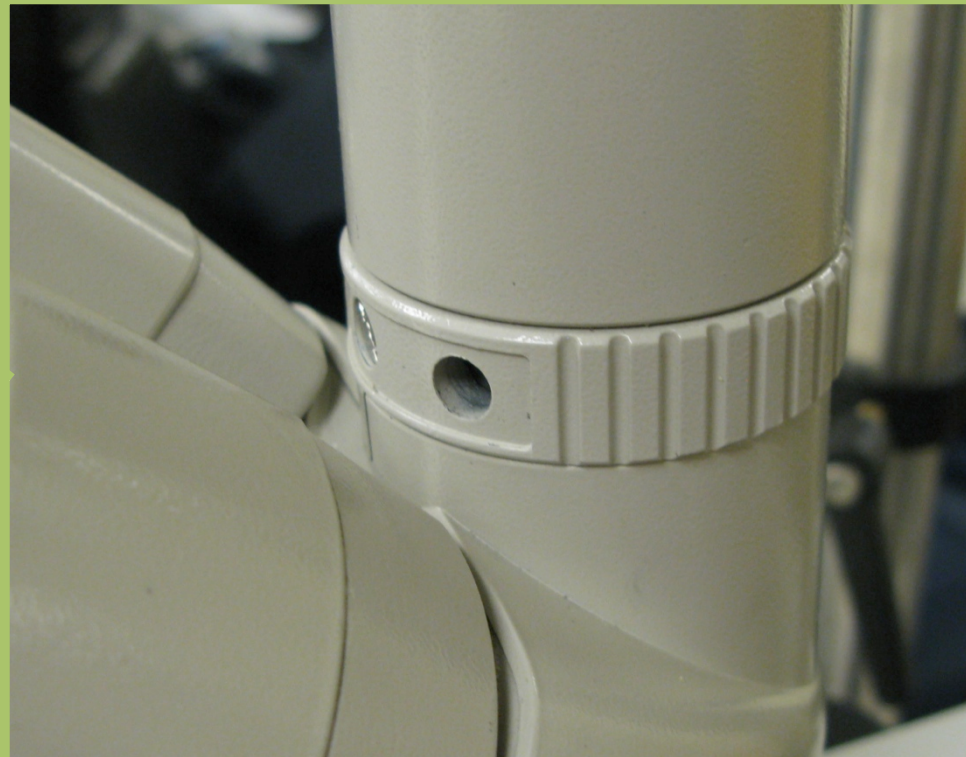

Course and fine focus can be found on either side of the microscope

Polarizing filter at base of light source rotate at the bottom of the unit

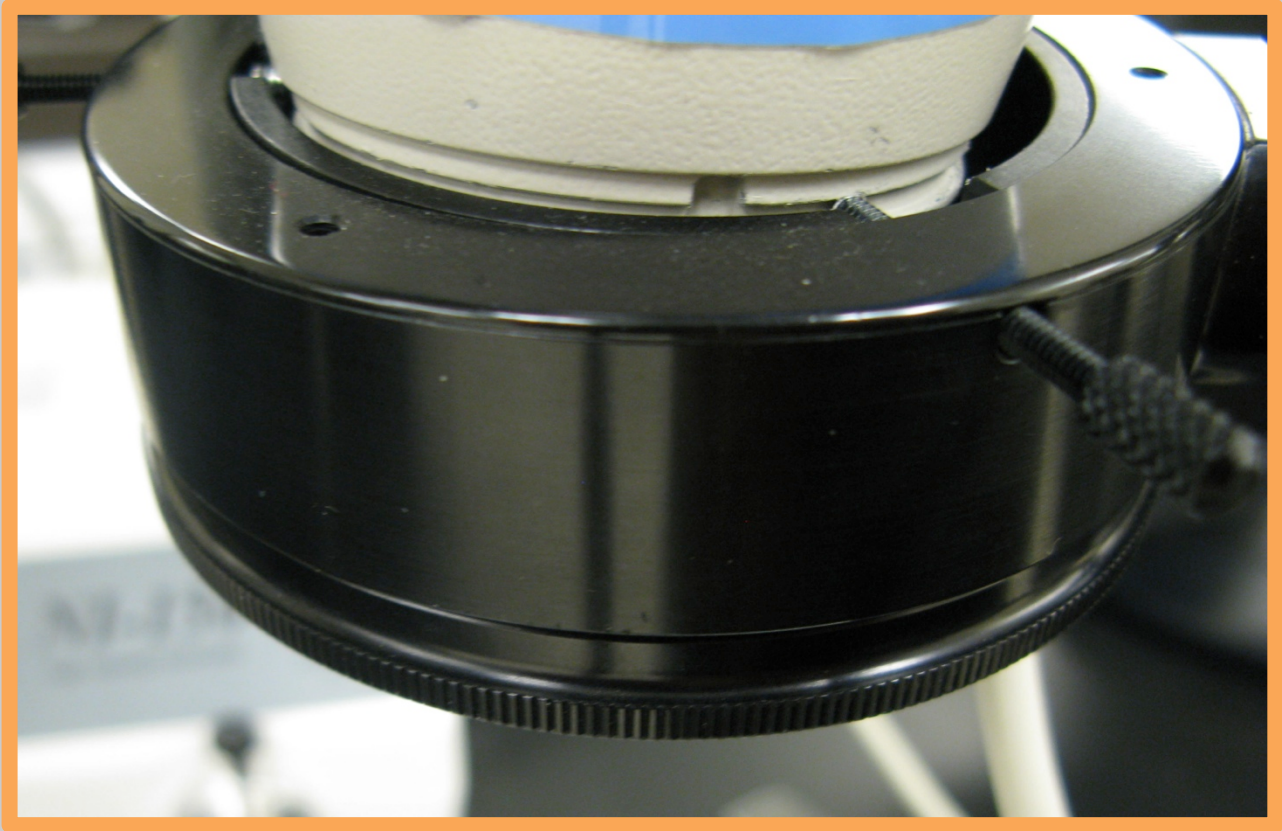

Power box for light source start at 0 power when turning on the unit